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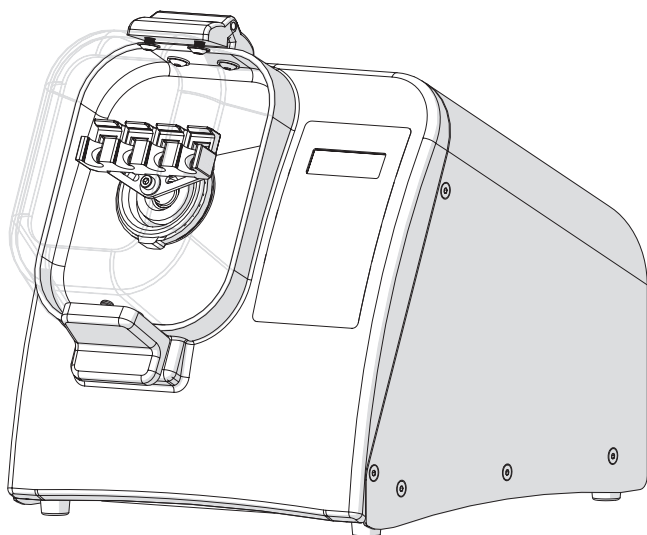
Homogeniser

Quick Start Guide




Catalog No	Description
432-0366	Mini Bead Mill 230V, 300W

Warning / Safety Information

- ⚠ **DO NOT** attempt to modify any part of this product.
- ⚠ **DO NOT** allow the machine to be submerged in any liquid.
- ⚠ **DO NOT** use in any setting other than an indoor laboratory.
- ⚠ **DO NOT** plug power cord into an incorrect outlet or subject it to an incorrect voltage.
- ⚠ Use this product only for its intended purpose.
- ⚠ **DO NOT** use attachments not recommended by the manufacturer.
- ⚠ **DO NOT** operate the product if it is damaged in any way.
- ⚠ Keep this product away from heated surfaces.
- ⚠ **DO NOT** modify the plug or cord that is provided. If the plug will not fit the outlet, have the proper outlet installed by a qualified electrician.
- ⚠ **DO NOT** operate the product with the safety ground disconnected.
- ⚠ **WARNING:** Reduce the risk of unintentional starting; make sure the speed switch is in the OFF position before plugging in the motor.
- ⚠ **WARNING:** Damaged or worn power cords should be repaired or replaced immediately by a qualified electrician.
- ⚠ **WARNING:** Improper connection of the equipment can result in a risk of electric shock.

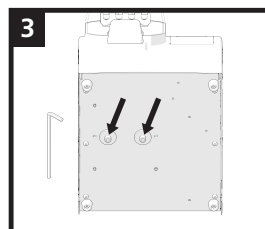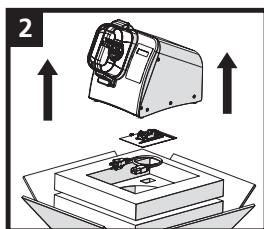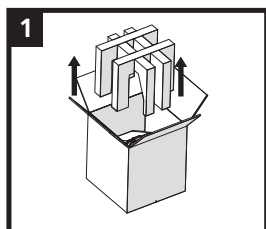
Unpacking


- ⚠ **WARNING:** Lift unit up and off of the box bottom.
DO NOT LIFT THE UNIT BY HOLDING THE LID.
Remove Tool Kit and Power Cord.

Remove the 2 screws from the bottom of the unit.

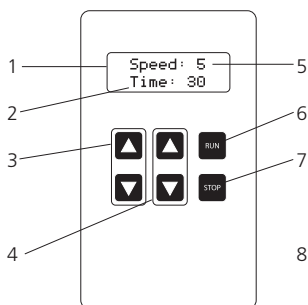General Operation

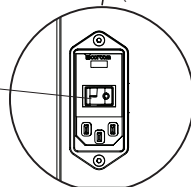1. Digital Display
2. Timer
3. Speed Control
4. Timer Control
5. Speed Setting
6. Run Button
7. Stop Button
8. On/Off switch

1. Power

- Plug in the power cord and move the On/Off switch on back of the unit to the On position.

2. Set Speed

- Adjust speed using up and down buttons.

3. Set Timer

- Adjust run time using up and down buttons from 1 sec. to 5 min.

4. Start

- Press Run button.



5. Pause/Stop

- Press Stop Button to Pause.

- Press Stop again to reset timer and cancel run

Speed Setting	Processing Power in G's	Processing Speed in m/s
1	30	1.5
2	60	2.5
3	90	3.7
4	120	4.4
5	150	5.0

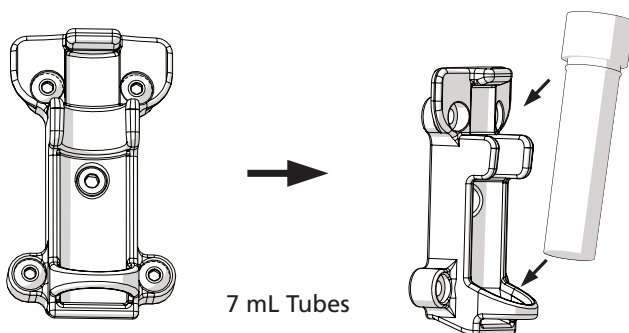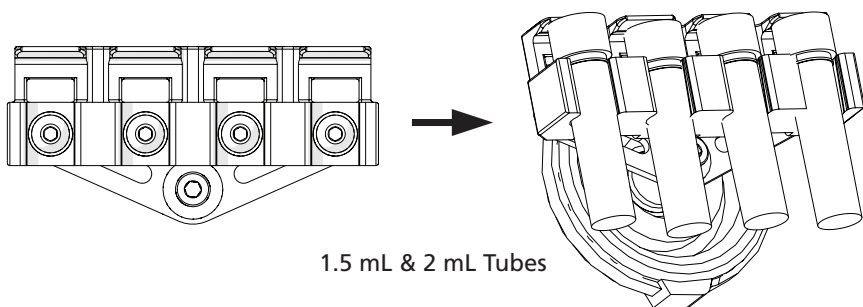Loading Sample Tubes

**For more details, please consult the full manual.
vwr.com**

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Simon László u. 4.
4034 Debrecen
Tel.: +36 (52) 521-130
Email: info.hu@vwr.com

India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

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Inde

VWR Lab Products Private Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road,
Peenya Post,
Bangalore, India - 560058
Tél. : +91-80-28078400
Courriel : vwr_india@vwr.com

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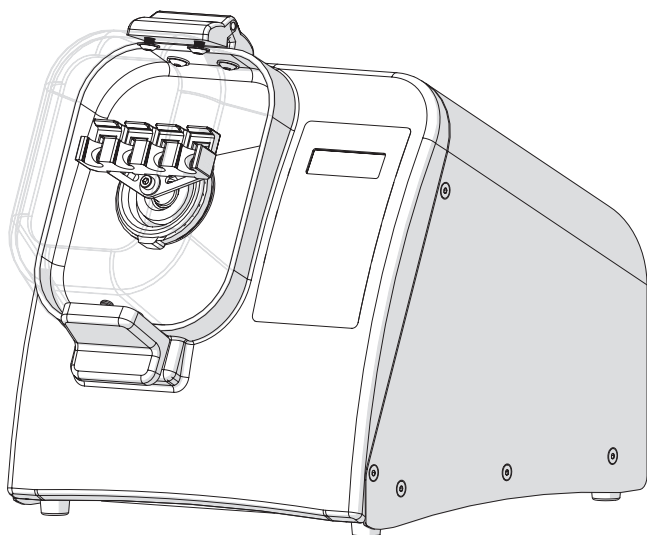
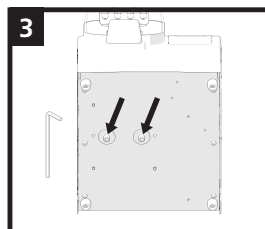
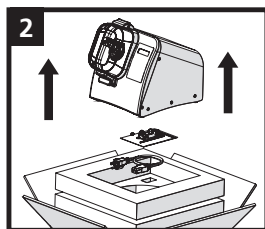
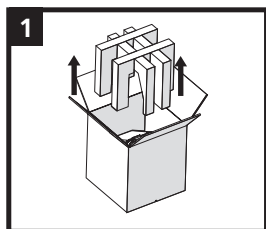
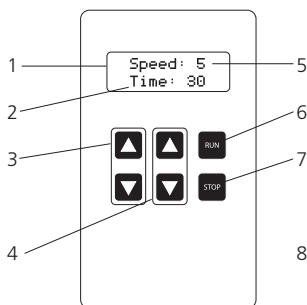
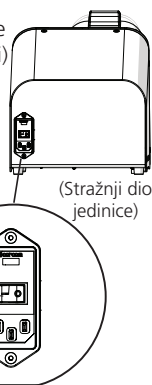
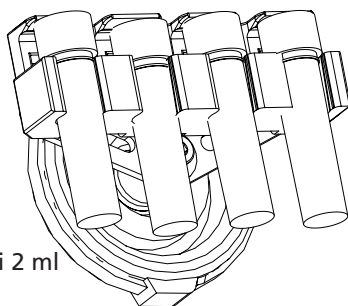
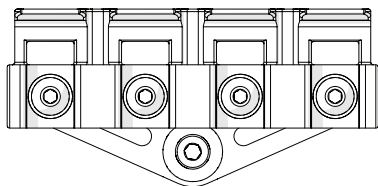
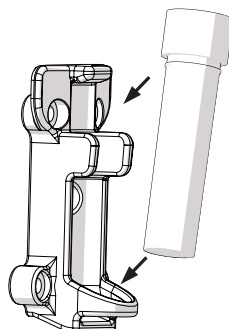
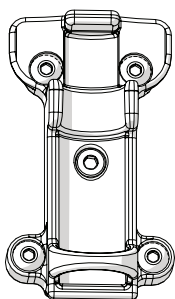
Indija

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road,
Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
E-mail: vwr_india@vwr.com

Irska / Sjeverna Irska

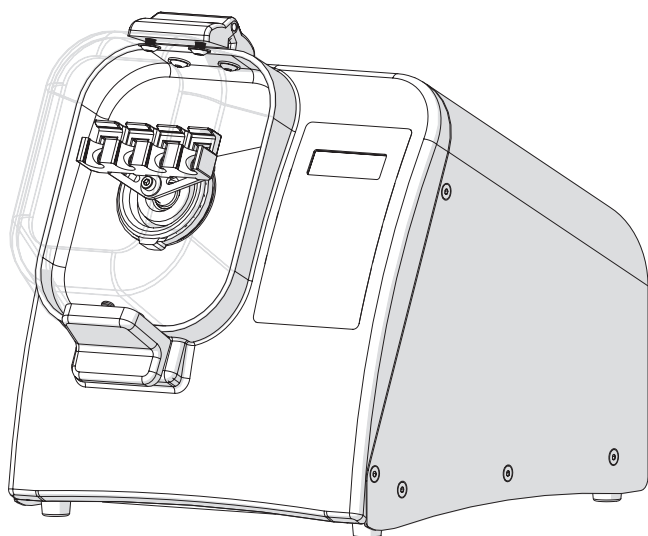
VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
E-mail sales.ie@vwr.com

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VWR International AB
VWR International GmbH