

Preparation of DNA Samples for Conventional DNA Electrophoresis

This protocol is for Preparation of DNA Samples for Conventional DNA Electrophoresis.

6x DNA Loading Dye, 6x Thermo Scientific MassRuler DNA Loading Dye (Cat #R0621), 6x Orange DNA Loading Dye (Cat #R0631), 6x Thermo Scientific TriTrack DNA Loading Dye (Cat #R1161) are all used according to below protocol:

1. Add 1 volume of 6x DNA loading dye to 5 volumes of DNA sample.
2. Mix well, spin down and load.

thermoscientific.com

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

North America
Technical Services:
ts.molbio@thermofisher.com
Customer Services:
cs.molbio@thermofisher.com
Tel 800 235 9880
Fax 800 292 6088

Europe and Asia
Technical Services:
ts.molbio.eu@thermofisher.com
Customer Services:
cs.molbio.eu@thermofisher.com
Tel 00800 2220888
Fax 00800 2220889

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific