

Magnetic Stand-96

Catalog Number AM10027

Pub. No. 4386455 Rev. D

Description	Quantity
Magnetic Stand-96	1

 **WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from thermofisher.com/support.

Product description

The Invitrogen™ Magnetic Stand-96 is designed for paramagnetic bead precipitation from standard 96-well U-bottom microplates and 0.2-mL PCR plates with no additional accessories. The stand has a white acetal copolymer base in a standard SBS 96-well microplate footprint. The use of 24 extremely strong permanent magnets guarantees easy and fast magnetic separation in as little as 30 seconds. The Magnetic Stand-96 can sit flat on a lab bench and in a microplate adapter on robotic workstations.

Using the Magnetic Stand-96

The Magnetic Stand-96 is ideal for use with the MagMAX™ *mirVana*™ Total RNA Isolation Kit, MagMAX™ DNA Multi-Sample Ultra Kit, and MagMAX™ Cell-Free DNA Isolation Kit. For more information on these kits, visit their web catalog pages at thermofisher.com.

IMPORTANT! We recommend cleaning the Magnetic Stand-96 with deionized water only; alcohol or alcohol-based cleaning agents may cause cracking of the stand material.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

The information in this guide is subject to change without notice.

DISCLAIMER

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. 4386455

Revision	Date	Description
D	27 April 2016	Updated the stand material and the list of compatible kits Format, style, and legal updates
C	27 June 2012	Baseline for this revision history

Important Licensing Information: This product may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Corporate entity: Life Technologies Corporation | Carlsbad, CA 92008 USA | Toll Free in USA 1 800 955 6288

©2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.