

# Reactive Oxygen Species (ROS) Fluorometric Assay Kit (Green)

Catalog Number EEA019 (96 tests)

Rev 4.0

For safety and biohazard guidelines, read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

## Product description

This kit can be used to measure reactive oxygen species (ROS) in fresh tissue and live cell samples.

Reactive oxygen species (ROS) are active chemical substances produced in the metabolic process of the body, including oxygen free radicals, hydrogen peroxide, and its downstream products, such as peroxides and hydroxides. ROS are both necessary and harmful to organisms, and are involved in cell growth, proliferation, development and differentiation, aging and apoptosis, as well as many physiological and pathological processes. Excessive ROS will lead to oxidative stress and oxidative damage of cells, further promoting the occurrence and development of many diseases, such as cancer, cardiovascular disease, and diabetes.

DCFH-DA (2,7-dichlorofluorescein diacetate) is a fluorescent probe without fluorescence that can freely cross the membrane. After entering the cell, it can be hydrolyzed by intracellular esterase to form DCFH (dichlorofluorescein). In the presence of reactive oxygen species (ROS), DCFH is oxidized to DCF (dichlorofluorescein) which is a strong green fluorescent substance that cannot penetrate the cell membrane. DCF has a maximum wave peak near the excitation wavelength of 502 nm and the emission wavelength of 525 nm, and the intensity is proportional to the level of intracellular reactive oxygen species.

## Contents and storage

Kit and components are shipped at -20°C protected from light. An unopened kit can be stored at -20°C for 12 months.

Components	Quantity (96 tests)
10 mmol/L DCFH-DA	0.1 mL
Positive Control	1 mL
Black Microplate	2 plates
Plate Sealer	4 pieces

## Required materials

- Double distilled water
- PBS (0.01 M, pH 7.4 )
- Serum-free medium
- Fluorescence Microplate reader (Ex/Em=488 nm/525 nm), Fluorescence Spectrophotometer (Ex/Em = 488 nm/525 nm), Flow Cytometry (Ex/Em = 488 nm/525 nm)
- Vortex mixer
- Micropipettor
- Water bath
- Incubator
- Centrifuge

## Procedural guidelines

---

**IMPORTANT!** Reagents are lot-specific. Do not mix or interchange different reagent lots from various kit lots.

---

Allow reagents to reach room temperature before use. Mix to redissolve any precipitated salts.

### Preparation of DCFH-DA working solution

Dilute the DCFH-DA with serum-free medium. The recommended working concentration is 0.1-20  $\mu\text{M}$ . (Note: DMSO is harmful to cells, so the dilution ratio must be more than 500.) Prepare the fresh solution before use.

### Preparation of positive control working solution

Dilute the positive control (contain 10 mM TBHP) with serum-free medium. The recommended working concentration of TBHP is 50-250  $\mu\text{M}$ . Prepare the fresh solution before use.

### The key points of the assay

- After the incubation of the probe, it is important to wash out residual probes that have not entered the cells, otherwise the background will be higher.
  - Avoid repeated freezing and thawing of DCFH-DA.
  - The time of detection is shortened as far as possible to reduce the experimental error.
- Set a positive control (positive control working solution) and a negative control (only cells without 10 mmol/L DCFH-DA working solution).

## Assay Protocol

### Detection of culture cell sample

#### Add the fluorescent probe

- a. Add DCFH-DA working solution to the cells. The DCFH-DA working concentration can be 0.1-20  $\mu\text{M}$  for different cells and treatment. A pre-experiment is suggested to determine the appropriate concentration. The total dilution ratio should be more than 500 in order to avoid effects of DMSO on cells. DMSO should be set as solution control.
- b. Incubate at 37°C for 30 min-few hours (generally 30~60 min). The incubation time is related to cell types, stimulation conditions, and DCFH-DA concentration.

**Cell collection:**

- a. Suspension cells: centrifuge the sample at 1000 g for 5-10 min and wash with serum-free medium for 2-3 times. Centrifuge and collect the cell precipitation for fluorescence detection.
- b. Adherent cells: digest the cells with 0.25% trypsin, then add medium that contains fetal bovine serum to terminate digestion, to prepare the cell suspension. Centrifuge at 1000 g for 5-10 min and collect cells, then wash with serum-free medium for 1-2 times. Centrifuge and collect cell precipitation for fluorescence detection.

**Fluorescence detection:**

- a. Re-suspend collected cells with serum-free medium for detection.
- b. Wavelength: the excitation wavelength is 488 nm, the emission wavelength is 525 nm. It can also be detected according to the fluorescence detection conditions of FITC.

**Notes:** the density of re-suspension cell is determined by cell fluorescence intensity. If fluorescence is strong (weak), then decrease (increase) the cell density, but cell density of all samples should be consistent.

**Detection of tissue sample****Preparation of single cell suspension****Method 1: using the single cell suspension instrument.****Method 2: enzyme digestion.**

- a. Take the tissue into pre-cooled PBS (0.01 M, pH 7.4) immediately and clean the blood and other contaminants. Remove the massive composition, fiber, fat, and blood vessels (except for specialized cells).
- b. Cut the tissue into about 1 mm<sup>3</sup> pieces with the ophthalmic scissors, then put these pieces to pre-cooled PBS (0.01 M, pH 7.4) to remove the cell debris.
- c. Add an appropriate amount of enzyme digestion, incubate in 37 °C water bath for 20~30 min, and gently oscillate the mixture intermittently.
- d. Stop the digestion with medium that contain fetal bovine serum. Filter the mixture to remove the tissue massive component with nylon mesh (300 mesh) and collect the cells. Centrifuge at 500 g for 10 min and discard the supernatant, then wash with PBS (0.01 M, pH 7.4) for 1~2 times. Re-suspend to prepare the single cell suspension solution. The cell amount should be no less than 10<sup>6</sup>.

### **Method 3: mechanical method.**

- a. The pretreatment is the same as step a and step b in the enzyme digestion method.
- b. Tight the nylon mesh (300 mesh) on a small beaker, then place the tissue pieces on the mesh and gently rub the tissue with ophthalmic scissor or erasing knife. Wash the tissue with PBS (0.01 M, pH 7.4) at the same time.
- c. Collect the cell suspension and centrifuge at 500 g for 10 min. Then discard the supernatant and wash with PBS (0.01 M, pH 7.4) for 1~2 times. Re-suspend to prepare the single cell suspension solution. The cell amount should be no less than  $10^6$

### **Add the fluorescent probe**

- a. Add DCFH-DA working solution to the cells. The DCFH-DA working concentration can be 0.1-20  $\mu\text{M}$  for different cells and treatment. A pre-experiment is suggested to determine the appropriate concentration. The total dilution ratio should be more than 500 in order to avoid effects of DMSO on cells. DMSO should be set as solution control.
- b. Incubate at 37°C for 30 min-few hours (generally 30~60 min). The incubation time is related to cell types, stimulation conditions, and DCFH-DA concentration.
- c. Collect the incubated single cell suspension, centrifuge at 1000 g for 5-10 min to collect cells. Wash with serum-free medium for 1~2 times. Centrifuge and collect the cell precipitation for fluorescence detection.

### **Fluorescence detection**

- a. Re-suspend collected cells with serum-free medium for detection.
- b. Wavelength: the excitation wavelength is 488 nm, the emission wavelength is 525 nm. It can also be detected according to the fluorescence detection conditions of FITC.

## **Notes**

- This kit is for research use only.
- Instructions should be followed strictly, changes of operation may result in unreliable results.
- The validity of kit is 12 months.
- Do not use components from different batches of kit.

- The density of re-suspension cell is determined by cell fluorescence intensity. If fluorescence is strong (weak), then decrease (increase) the cell density, but cell density of all samples should be consistent.
- Fluorescent substances are sensitive to light and should avoid light during detection.
- DCF is easy to be quenched, and the samples after incubation must be detected within 2 hours.
- Results were expressed as fluorescence intensity or geometric average fluorescence intensity (flow cytometry).
- When using flow cytometry, in order to avoid the interference of cell debris and dead cells on the experimental results, it is necessary to eliminate them.
- Set a positive control (positive control working solution) and a negative control (only cells without 10 mmol/L DCFH-DA working solution).
- The timing of adding DCFH-DA or incubation time depends on whether the intracellular reactive oxygen species can be detected successfully. DCFH-DA can be added in advance or at the same time if the drug treatment time is short ( $< 2$  h) or the predicted ROS is weak. Conversely, DCFH-DA can be added later if the drug treatment time is long ( $> 6$  h) or predicted ROS is strong.

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at

**[www.thermofisher.com/us/en/home/global/terms-and-conditions.html](http://www.thermofisher.com/us/en/home/global/terms-and-conditions.html)**. If you have any questions, please contact Life Technologies at

**[www.thermofisher.com/support](http://www.thermofisher.com/support)**.

Corporate entity: Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

The information in this guide is subject to change without notice.

### DISCLAIMER

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

© 2025 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.