CAMP TEST- S

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1. Christie R., Atkins N.E., Munch- Peterson E. "A note on a lytic phenomenon shown by group B streptococci". Aug. J. Biol. 22: 197-200 (1994).
2. Brzin B. and H.P.R. Seeliger (1975). "A brief note on the CAMP phenomenon in listeria". In problems of Listeriosis (M. Woodbine, Ed), Leichestor University press.
3. ISO 11290-1:1196. Microbiology of food and animal feeding stuffs – Horizontal method for the detection and enumeration of *Listeria monocytogenes* – Part 1: Detection method.

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CAMP TEST- S

Rapid test for identification of *Listeria* spp.

DESCRIPTION

CAMP TEST- S is constituted by paper discs impregnated with *Staphylococcus aureus* ATCC 25923 used to estimate haemolytic power of group B streptococci and listeriae.

PRINCIPLE

CAMP factor is an extracellular substance produced by group B streptococci and some strains of *Listeria* which increases lysis of red cells by beta- lysin of *Staphylococcus aureus*. *Staphylococcus aureus* used for the impregnation is a viable and stabilized microorganism resulting from the collection of the American Type of Culture Collection (ATCC).

TECHNIQUE

1. Take a disc from a tube and put it down on the surface of a plate of Tryptic Soy Agar + 5% Sheep Blood (ref. 11037).
2. Moisten the disc with a drop of sterile physiological solution.
3. Touch with a sterile loop the bacterial colony to examine, mark out one or two lines near the disc, caring not to touch the edge of the disc.
4. Incubate the plate at 36+/-1°C for 24 ore.

INTERPRETATION OF RESULTS

Observe for the presence of a haemolytic zone near the disc and interpret according to the scheme:

ENHANCED ZONE OF HAEMOLYSIS NEAR THE DISC	POSITIVE TEST
NO HAEMOLYSIS NEAR THE DISC	NEGATIVE TEST

Scheme of CAMP TEST reaction for *Listeria* spp:

MICROORGANISM	BETA-HAEMOLYSIS	CAMP TEST- S
<i>Listeria monocytogenes</i>	+	+
<i>Listeria innocua</i>	-	-
<i>Listeria ivanovii</i>	+	-
<i>Listeria seelegeri</i>	+	+
<i>Listeria welshimeri</i>	-	-
<i>Listeria grayi</i> subspecies <i>grayi</i>	-	-
<i>Listeria grayi</i> subspecies <i>murrayi</i>	-	-
<i>Jonesia denitrificans</i>	-	-

STORAGE AND TRASPORT CONDITIONS

2-8°C away from light, until the expiry date on the label. However, our stability studies have shown that the storage or transport at 18-25°C for 4 days, or at 35-39°C for 48 hours, do not alter in any way the performance of the product. Eliminate if signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is designed for *In vitro* diagnostic use and must be used only by properly trained operators.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to the national and local regulations in force.

REFERENCES

1. Christie R., Atkins N.E., Munch- Peterson E. "A note on a lytic phenomenon shown by group B streptococci". *Aug. J. Biol.* 22: 197-200 (1994).
2. Brzin B. and H.P.R. Seeliger (1975). "A brief note on the CAMP phenomenon in listeria". In *problems of Listeriosis* (M. Woodbine, Ed), Leichestre University press.
3. ISO 11290-1:1196. Microbiology of food and animal feeding stuffs – Horizontal method for the detection and enumeration of *Listeria monocytogenes* – Part 1: Detection method.

LIOFILCHEM® S.r.l.

Via Scozia, Zona Ind.le - 64026, Roseto degli Abruzzi (TE) - ITALY
Tel +39 0858930745 Fax +39 0858930330 Website: www.liofilchem.net E-mail: liofilchem@lioilchem.net

PRODUCT SPECIFICATIONS

NAME

CAMP TEST- S

PRESENTATION

30 paper disc in cartridge

STORAGE

2-8°C

PACKAGING

Ref.	Content	Packaging
88021	30 discs	<ul style="list-style-type: none"> • 1 cartridge in a plastic bottle • 1 plastic bottle in a cardboard box

USE

CAMP TEST- S is constituted by paper discs impregnated with *Staphylococcus aureus* ATCC 25923 used to estimate haemolytic power of group B streptococci and listeriae

TECHNIQUE

Refer to technical sheet of the product

SHELF LIFE

1 year

QUALITY CONTROL

1. Control of general characteristics, label and print
2. Microbiological control
Incubation Conditions: 18-24 h at 35 ± 2°C

Microorganisms		Enhanced Zone of Haemolysis
<i>Listeria monocytogenes</i>	CIP 7836	+
<i>Listeria innocua</i>	CIP 8011	-
<i>Listeria welshimeri</i>	CIP 8149	-

TABLE OF SYMBOLS

LOT	Batch code	IVD	<i>In vitro</i> Diagnostic Medical Device		Manufacturer		Use by		Fragile, handle with care
REF	Catalogue number		Temperature limitation		Contains sufficient for <n> tests		Caution, consult instruction for use		Keep away from heat sources

LIOFILCHEM® S.r.l.

Via Scozia, Zona Ind.le - 64026, Roseto degli Abruzzi (TE) - ITALY
Tel +39 0858930745 Fax +39 0858930330 Website: www.liofilchem.net E-mail: liofilchem@liofilchem.net