


Spin Columns and Elution Tubes

Catalog Number AM10065

Pub. No. 4386458 Rev. C

Contents	Quantity	Storage conditions
Spin Columns and 2 mL Elution Tubes	100 each	Ambient

 **WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from www.lifetechnologies.com/support.

Product description

Spin Columns and Elution Tubes can be used as a simple way to recover DNA from agarose gels after electrophoresis, during plasmid purification procedures, and in poly(A) selection of mRNA.

These Spin Columns and Elution Tubes may also be used to collect Oligo(dT) Cellulose (Cat. no. AM10020) bound with mRNA.

Recovery of DNA from agarose gels

1. Check that the Spin Column/Elution Tube Assembly fits into your microcentrifuge when spun.
2. Excise the DNA or RNA band of interest from the gel.
3. Place the gel slice in a spin column, suspended in the provided 2 mL Elution Tubes, and centrifuge for 5 min at top speed. The nucleic acid will be forced through the column and into the tube below while the agarose is left behind. The nucleic acid is then ready for manipulations, such as restriction enzyme digestion, ligation, or reverse transcription.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

For Research Use Only. Not for use in diagnostic procedures.

The information in this guide is subject to change without notice.

DISCLAIMER: LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

NOTICE TO PURCHASER: LIMITED USE LABEL LICENSE: Research Use Only : The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

© 2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

For support visit lifetechnologies.com/support or email techsupport@lifetech.com

lifetechnologies.com

18 October 2012

