

Performance characteristics

Standard curve (example)

The following data were obtained for the various standards over the range of 0-92 U/mL SOD.

Standard SOD (U/mL)	Optical Density (450 nm)
92.0	0.124
46.0	0.214
23.0	0.355
11.5	0.500
5.8	0.652
2.9	0.774
1.4	0.860
0.0	0.939

Intra-assay precision

Three samples were assayed in replicates of 20 to determine precision within an assay.

Parameters	Sample 1	Sample 2	Sample 3
Mean (U/mL)	25.5	15.5	7.4
%CV	2.0	3.0	1.4

CV = Coefficient of Variation

Inter-assay precision

Three samples were assayed 20 times in duplicate by five operators to determine precision between assays.

Parameters	Sample 1	Sample 2	Sample 3
Mean (U/mL)	23.7	13.5	6.6
%CV	8.3	10.9	15.3

CV = Coefficient of Variation

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Catalog Number

Batch code

Temperature limitation

Use by

Manufacturer

Consult instructions for use

Caution, consult accompanying documents

Manufacturer: Life Technologies Corporation | 7335 Executive Way | Frederick, MD 21704 | USA

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2020 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

Superoxide Dismutase (SOD) Colorimetric Activity Kit

Catalog Number EIASODC (192 tests)

Pub. No. MAN0019052 Rev A.0

For safety and biohazard guidelines, see the “Safety” appendix in the *ELISA Technical Guide* (Pub. no. MAN0006706). Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

Product description

The Superoxide Dismutase (SOD) Colorimetric Activity Kit is designed to measure all types of SOD activity (i.e., Cu/Zn, Mn, and Fe superoxide dismutases) in a variety of samples.

This assay measures the activity of SOD in serum, plasma, tissue lysates, cell lysates, and erythrocyte lysates. The assay was characterized with human SOD, but is expected to measure SOD activity in samples from other species.

Reactive oxygen species are produced during normal aerobic metabolism. These free radicals are removed or converted to other products *in vivo* by the activity of specific enzymes including a variety of superoxide dismutases which play important roles in scavenging O₂⁻. Cu/Zn SOD (SOD1) is the major intracellular superoxide dismutase. MnSOD (SOD2) is the mitochondrial superoxide dismutase, and has similar properties to FeSOD found in some plants and bacteria. EC-SOD (SOD3) is the primary extracellular superoxide dismutase, and can be detected in serum, plasma, ascites, or synovial fluids.

Contents and storage

Kit and components are shipped at –20°C. Upon receipt, store the kit at –20°C. Once open, store the kit at 4°C and use within 2 weeks. Store the Superoxide Dismutase Standard at –20°C or lower after opening.

Components	Quantity
Superoxide Dismutase Standard; 9,200 U/mL bovine superoxide dismutase	50 µL
Assay Buffer; containing detergents, stabilizers, and dye	50 mL
Clear 96-well Half Area Plate	2 plates
Substrate Concentrate (10X)	1.1 mL
Substrate Diluent; keep tightly capped	12 mL
Xanthine Oxidase Buffer; containing detergents and stabilizers	6 mL
Xanthine Oxidase Concentrate (25X)	225 µL

Materials required but not supplied

- Microtiter plate reader with software capable of measurement at or near 450 nm
- Calibrated adjustable precision pipettes and glass or plastic tubes for diluting solution
- (*Optional*) Potassium cyanide stock solution

Procedural guidelines

Reagents are lot-specific. Do not mix or interchange different reagent lots from various kit lots.

Sample preparation guidelines

- Collect samples in pyrogen/endotoxin-free tubes.
- Freeze samples after collection if samples will not be tested immediately. Avoid multiple freeze-thaw cycles of frozen samples. Thaw completely and mix well (do not vortex) prior to analysis.
- Avoid the use of hemolyzed or lipemic sera.
- If large amounts of particulate matter are present in the sample, centrifuge or filter sample prior to analysis.

Keep samples on ice to maintain enzyme activity.

Prepare samples

Sample	Procedure
Serum/ plasma	<p>Collect samples</p> <ul style="list-style-type: none">• Collect blood in serum tubes. Centrifuge at 700 × <i>g</i> for 15 minutes at 4°C. Aspirate the supernatant. Assay immediately or freeze at –80°C.• Collect plasma in heparin or EDTA tubes. Centrifuge at 700 × <i>g</i> for 15 minutes at 4°C. Aspirate the pale yellow supernatant. Assay immediately or freeze at –80°C. <p>NOTE: Some serum and plasma samples may contain significant hemoglobin concentrations and may require blanking prior to adding 1X Xanthine Oxidase (see Assay procedure, page 3).</p>
Erythrocytes (RBCs)	<ol style="list-style-type: none">1. Eyrthrocytes can be lysed by taking the pelleted RBCs from plasma as described in the previous section and adding 4 volumes of ice cold deionized water.2. Centrifuge at 10,000 × <i>g</i> for 15 minutes at 4°C to remove debris. <p>NOTE: Lysed RBCs will exhibit high background color and require blanking prior to adding 1X Xanthine Oxidase (see Assay procedure, page 3).</p>

Prepare samples, continued

Sample	Procedure
Cell lysate	<div><div>1.</div><div>Prepare cell pellet.<div><div>•</div><div>For cells in suspension, centrifuge >1 × 10⁶ cells at 250 × <i>g</i> for 10 minutes at 4°C, then discard the supernatant.</div><div>•</div><div>For adherent cells, rinse the culture dish with PBS, then harvest >1 × 10⁶ cells by gentle trypsinization. Transfer to a tube on ice and centrifuge at 250 × <i>g</i> for 10 minutes at 4°C and discard the supernatant.</div></div></div><div><div>2.</div><div>Resuspend the cell pellet in ice-cold PBS and transfer to a microcentrifuge tube on ice. Centrifuge the cells at 250 × <i>g</i> for 10 minutes at 4°C, discard the supernatant, and place the tube on ice.</div></div><div><div>3.</div><div>Homogenize or sonicate the pellet in 0.5–1 mL of PBS per 100 mg of cells. Centrifuge at 1,500 × <i>g</i> for 10 minutes at 4°C.</div></div><div><div>4.</div><div>Collect the supernatant and assay immediately, or store at -80°C. To measure cytosolic (SOD1, Cu/Zn) and/or mitochondrial SOD (SOD2, Mn) centrifuge supernatants at 10,000 × <i>g</i> for 15 minutes at 4°C. The supernatants will contain the cytosolic SOD and the cell pellets will contain mitochondrial SOD. To determine Mn SOD (SOD2) activity treat samples with potassium cyanide to a final concentration of 2 mM. Addition of cyanide will inactivate other SOD enzymes</div></div></div>
Tissue lysate	<div><div>1.</div><div>Wash tissue thoroughly with ice cold PBS prior to processing.</div></div> <div><div>2.</div><div>Homogenize tissue as described in previous section for cell lysate, steps 3 and 4.</div></div>

Dilute samples


Sample concentrations should be within the range of the standard curve. Because conditions may vary, each investigator should determine the optimal dilution for each application.

- Dilute **cell and tissue lysate** samples ≥1:4 in Assay Buffer.
- Dilute **serum and plasma** samples ≥1:5 in 1X Assay Buffer.
- Dilute **erythrocyte** samples ≥1:100 in Assay Buffer.
- Use all samples within **2 hours** of dilution.

Dilute standards

Note: Use glass or plastic tubes for diluting standards.
Note: One unit of SOD is the amount of enzyme causing half the maximum inhibition of the oxidation of 7.5 mM NADH in the presence of EDTA, manganese ions, and mercaptoethanol at 23°C and pH 7.4 over 15 minutes.

1. Add 10 µL SOD Standard to one tube containing 990 µL Assay Buffer and label as 92 U/mL SOD.
2. Add 100 µL Assay Buffer to each of 7 additional tubes labeled as follows: 46, 23, 11.5, 5.8, 2.9, 1.4, and 0 U/mL SOD.
3. Make serial dilutions of the standard as described below in the dilution diagram. Mix thoroughly between steps.
4. **Use the standards within 2 hours of preparation.** Store remaining reconstituted standard in frozen aliquots at –20°C or lower.

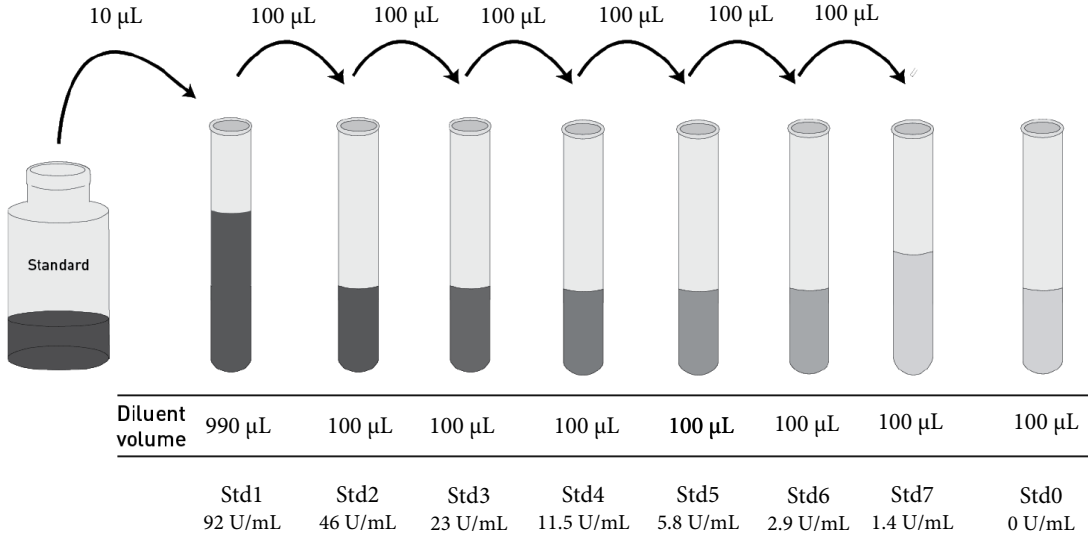

Prepare 1X Xanthine Oxidase

Vortex Xanthine Oxidase Concentrate prior to pipetting. Pipet from the bottom of the tube.

Reagent	½ plate	1 plate	1½ plates	2 plates
Xanthine Oxidase	50 µL	100 µL	150 µL	200 µL
Xanthine Oxidase Buffer	1.2 mL	2.4 mL	3.6 mL	4.8 mL
Total volume	1.25 mL	2.5 mL	3.75 mL	5 mL

Prepare 1X Substrate

Vortex the vial of Substrate Concentrate prior to pipetting.

Note: Keep Substrate Diluent and 1X Substrate tightly capped when not in use.

Reagent	½ plate	1 plate	1½ plates	2 plates
Substrate Concentrate	250 µL	500 µL	750 µL	1 mL
Substrate Diluent	2.25 mL	4.5 mL	6.75 mL	9 mL
Total volume	2.5 mL	5 mL	7.5 mL	10 mL

Assay procedure


Allow all reagents to reach room temperature before use. Mix all liquid reagents prior to use. **Total assay time is 20 minutes.**

IMPORTANT! Perform a standard curve with each assay.



Add sample and substrate

- a. Add 10 µL of standards or diluted samples (see page 2) to the appropriate wells.
- b. Add 50 µL of 1X Substrate into each well.

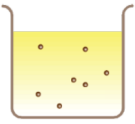

Add chromogenic detection reagent ^[1]

- a. Add 25 µL of 1X Xanthine Oxidase into each well.
- b. Incubate for 20 minutes at room temperature.

[1] Bright yellow colored samples can interfere with the high sensitivity format assay, and may require blanking prior to adding 1X Xanthine Oxidase.

To blank the plate, read the absorbance at 450 nm before adding the chromogenic detection reagent. Subtract the absorbance values from the blanking step from the corresponding absorbance measured for each sample.

Read the plate and generate the standard curve