

Self-circularization of Linear DNA

This protocol is for the Self-circularization of Linear DNA

1. Prepare the following reaction mixture:

Linear DNA	10-50 ng
10X T4 DNA Ligase buffer	5 μ l
T4 DNA Ligase	5 u
Water, nuclease-free	to 50 μ l
Total volume	50 μl

2. Mix thoroughly, spin briefly and incubate:
 - sticky-ends for 10 min at 20°C,
 - blunt-ends for 1 hour at 22°C.
3. Use up to 5 μ l of the mixture for transformation of 50 μ l chemically competent cells and 1-2 μ l per 50 μ l of electrocompetent cells.

Note

- The electrotransformation efficiency may be improved by:
 - heat inactivation of T4 DNA Ligase at 65°C for 10 min or at 70°C for 5 min,
 - purification of DNA, using GeneJET™ PCR Purification Kit (#K0701), or by chloroform extraction.
- The overall number of transformants may be increased by extending the reaction time to 1 hour.

IMPORTANT NOTES

- Polyethylene glycol (PEG) greatly increases the ligation efficiency of blunt-end DNA ligation. The recommended concentration of PEG 4000 in the ligation reaction mixture is 5% (w/v).
- Do not exceed the recommended amount of T4 DNA Ligase in the reaction mixture.
- Binding of T4 DNA Ligase to DNA may result in a band shift in agarose gels. To avoid this, incubate samples with 6X Loading Dye & SDS Solution (#R1151) at 70°C for 5 min and chill on ice prior to loading.
- For efficient transformation, the volume of the ligation reaction mixture should not exceed 10% of the competent cell volume.

thermoscientific.com

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

North America
Technical Services:
 techservice.genomics@thermofisher.com
Customer Services:
 customerservice.genomics@thermofisher.com
 Tel 800 235 9880
 Fax 800 292 6088

Europe and Asia
Technical Services:
 techservice.emea.genomics@thermofisher.com
Customer Services:
 customerservice.emea.genomics@thermofisher.com

Thermo
 S C I E N T I F I C

Part of Thermo Fisher Scientific